

# Psychtoolbox-3 (PTB3)

## Simple Display Calibration for Grayscale Stimuli

### Using a Minolta CS100A Photometer

(Frank Schieber, University of South Dakota, [schieber@usd.edu](mailto:schieber@usd.edu) June 13, 2007)

#### Abstract

This document briefly describes a set of PTB3/MATLAB scripts developed to perform simple grayscale calibration and linearization of an RGB monitor. See the built-in help and comments in each script for additional information. This code was developed and tested in a Windows environment. Mac users should be able to implement the code with little or no modification assuming that the serial I/O routines function properly (i.e., `openCS100A.m`, `readCS100A.m` and `closeCS100A.m`).

#### Development and Test Environment

OS: Windows XP Pro (SP2)  
MATLAB: 7.0.4.365 (R14) Service Pack 2  
Graphics card: ATI Radeon 9600  
Photometer: Minolta CS100A (Requires special cable to connect to RS232 serial port)  
Modify `openCS100A/readCS100A/closeCS100A` if you need to substitute another photometer. Only luminance readings are used here.

#### Files needed (over and above PTB3 distribution)

<u>Filename</u>	<u>Description</u>
<code>openCS100A.m</code>	open RS232 serial port connection to CS100A meter
<code>readCS100A.m</code>	read luminance and chromaticity from CS100A meter
<code>closeCS100A.m</code>	close persistent connection to CS100A meter
<code>phase1.m</code>	collects luminance measurements from display using a <u>linear gamma table</u> and plots results [results saved in <i>phase1_photometry.mat</i> file]
<code>phase2.m</code>	generates best-fit power function to the luminance data collected by <code>phase1.m</code> and then builds an inverse gamma table to compensate for grayscale display nonlinearities. Generates summary plots and saves results in <i>phase2_photometry.mat</i>

<u>Filename</u>	<u>Description</u>
phase3.m	installs inverse gamma table generated by phase2.m and measures screen luminance at every possible grayscale value [0:255]. Results are saved in <i>phase3_photometry.mat</i> .
phase4.m	generates a graphical summary of all results from calibration phases 1-3. Check here to see how well the inverse gamma table succeeded in “linearizing” the grayscale display output

### Calibration procedure in a nutshell

#### Step 0.

The first thing that you need to do is make sure that the photometer and MATLAB can successfully communicate with one another. The Minolta CS100A photometer requires a special communications cable which plugs into the photometer at one end and a standard RS232 cable at the other end. The RS232 (serial) cable is then connected to the COM1 communications port on the computer. To coerce the CS100A to enter RS232 COMMUNICATIONS mode, you need to hold down the “F” key on the photometer while you transition the power switch from the “off” to “on” positions (A small letter “c” will appear on the LCD display to verify COMM mode). Next, execute the **openCS100A** script to establish a serial port connection between MATLAB and the photometer (The script returns a zero if successful). Once this connection is established, you can read the photometer as follows:


[luminance, x, y, status] = **readCS100A**

When finished using the photometer, good programming practice mandates that you close the serial port connection using the **closeCS100A** script.

#### Step 1.

Now that you have verified a working connection to the photometer you can begin taking some preliminary measurements to characterize the photometric properties of your display system. Setup the photometer on a tripod and focus it on the display screen. Execute the **phase1.m** script. This script opens a PTB3 screen and prompts you to press the Enter key on the keyboard to begin. However, first you should take the opportunity to aim the photometer on the fixation point (“+”) near the center of the display. When you press the enter key, a 10 sec (default) countdown will begin to give you time to turn out the lights (and exit the room). At this point, the script generates and uploads a linear gamma table to the graphics card and then sequentially increments the luminance index of the calibration target from 0 to 255 (in steps of size 8) while the photometer collects

luminance readings under program control. Feedback regarding the current index is provided in the lower left corner of the display. When the procedure is completed, a luminance-by-index curve is plotted and saved to *phase1\_photometry.mat* (see Figure 1).

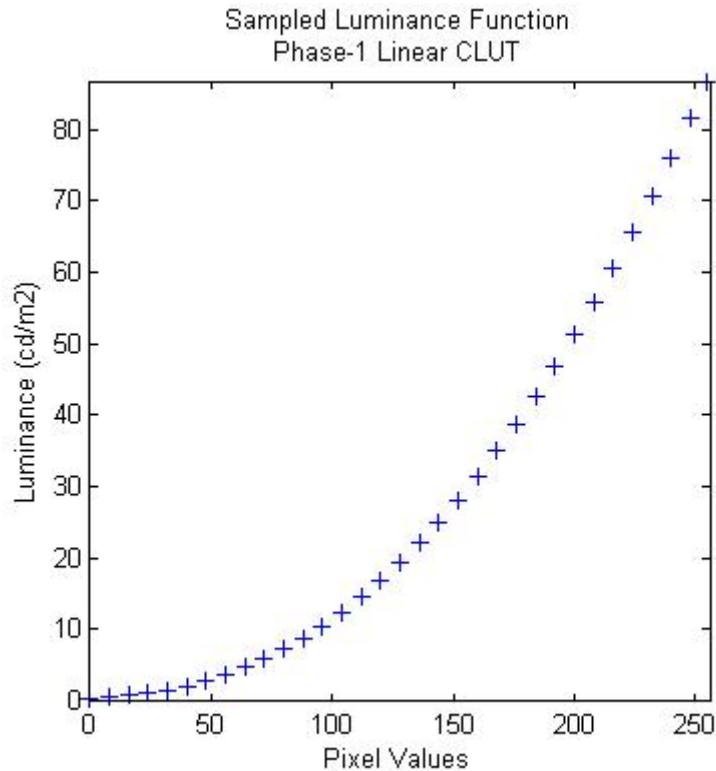


**Figure 1.** Sample output from phase1.m script.


### Step 2.

Now that you have characterized the nonlinearity that exists between a linear gamma table in your display adapter and the actual luminance output, you can execute the **phase2.m** script to generate a best fitting power function to the observed luminance values and then use the resulting parameters to build an inverse gamma table<sup>1</sup> to “linearize” the luminance output of your display. The results of this process are summarized in a MATLAB plot and saved to *phase2\_photometry.mat* (See Figures 2 and 3).

<sup>1</sup> As currently implemented, the phase2.m script does not double-check the inverse gamma table for violations of monotonicity.

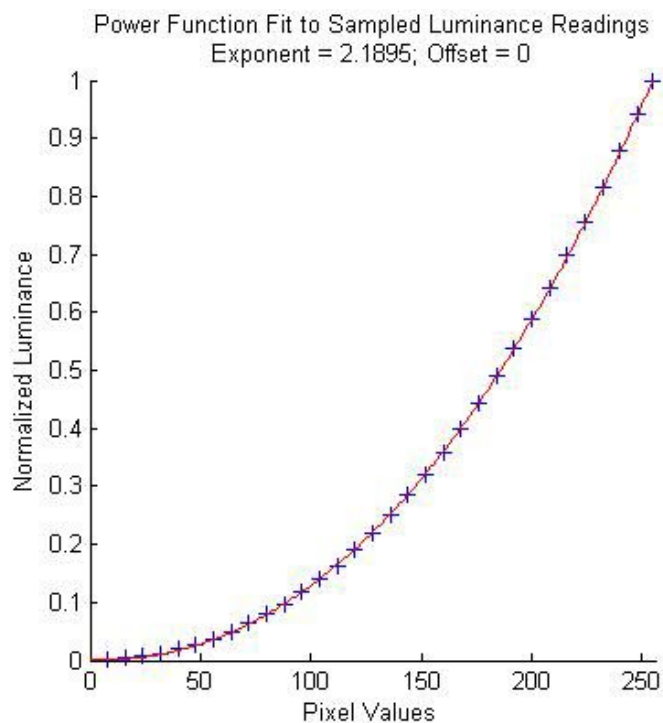


**Figure 2.** Best-fit power function to luminance data collected with linear gamma table.

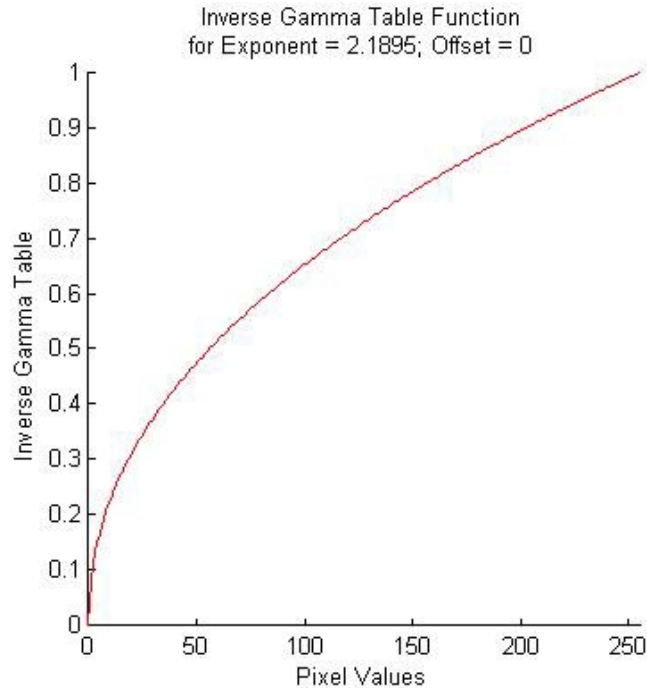

**Figure 3.** Inverse gamma table generated by phase2.m to compensate for non-linearity depicted in Figures 1 and 2.

### Step 3.

Now that you have generated and saved an inverse gamma table designed to compensate for the luminance non-linearity revealed in steps 1-2, it's time to load the inverse gamma table and evaluate how well it achieves its purpose. Execute **phase3.m** to load the inverse gamma table and collect luminance data from your display. This script has the same “look and feel” as *phase1.m* (described above) except that it takes much longer to run since it samples all possible index values from 0 thru 255. The results are automatically saved to *phase3\_photometry.mat* and can be plotted by executing the **phase4.m** script. Figure 4 depicts an example of the output generated by *phase4.m* and reveals that the inverse gamma table did a reasonably good job of “linearizing” the luminance output of the display subsystem (see the long URL at the bottom of this page for Jenny Read’s “brute force” recursive technique for achieving optimal gamma table correction results).

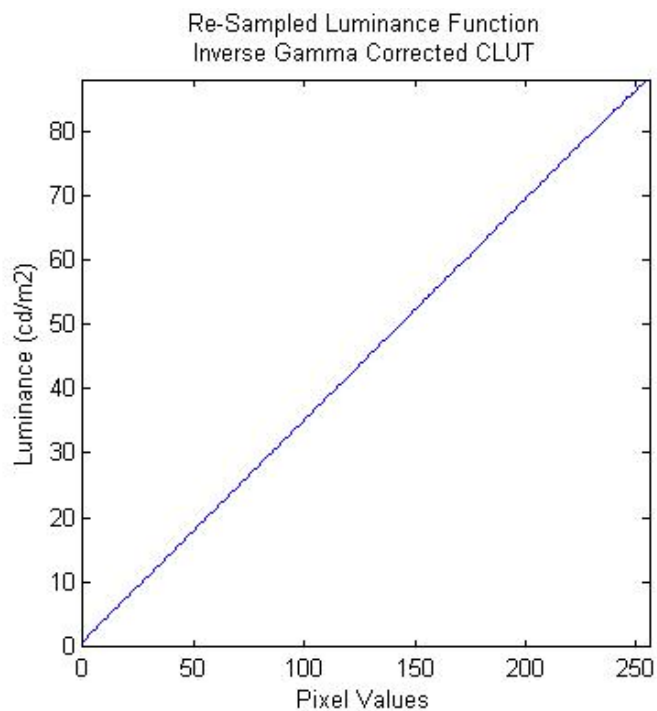

**Figure 4.** Linearized luminance output achieved by inverse gamma table correction.

Jenny Read's Display Linearization Web Page (Highly recommended)

<http://www.staff.ncl.ac.uk/j.c.a.read/index.php?location=research&sub=labsetup&file=research/labsetup/gammacorrection.html>